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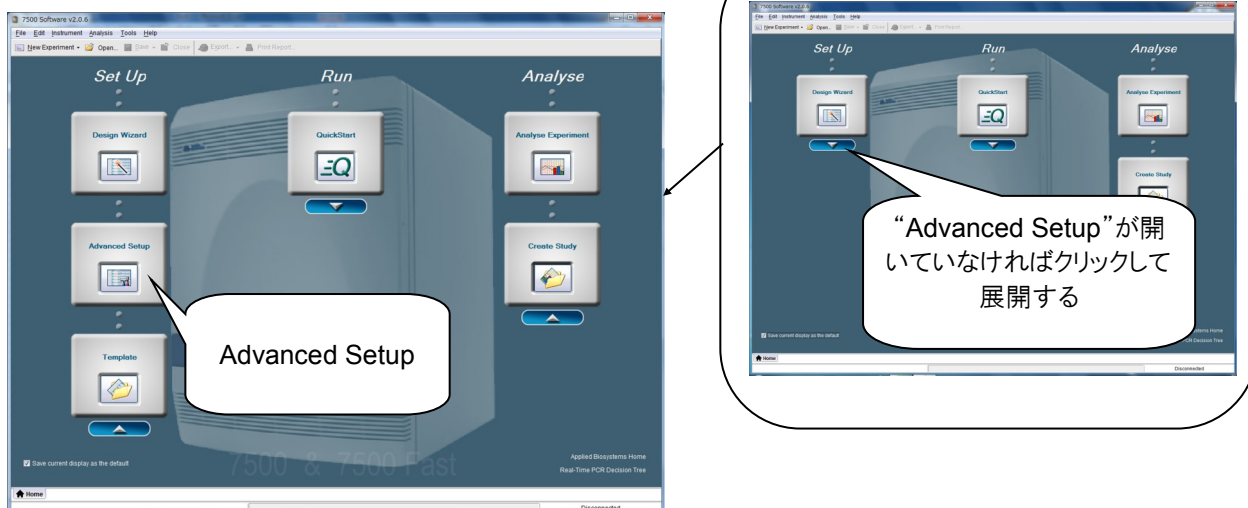
## 2) “Advanced Setup” を開く

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The screenshot shows the 'Experimental Properties' window in 7500 Software v2.0.6. The window title is 'Experiment: 20151214name Type: Standard Curve Reagents: SYBR® Green Reagents'. The 'Experiment Properties' section includes fields for 'Experiment Name' (20151214name), 'Barcode (Optional)', 'User Name (Optional)', and 'Comments (Optional)'. A callout points to the 'Experiment Name' field with the text: 'Date + free word Ex) 20151224ocu'. The 'Which instrument are you using to run the experiment?' section has a dropdown menu set to '7500 Fast (96 Wells)'. A callout points to this dropdown: '装置の選択です。必ず 7500Fastを選択してください。'. The 'Which reagent do you want to use to detect the target sequence?' section has a dropdown menu set to 'SYBR® Green Reagents'. A callout points to this dropdown: '使用している蛍光試薬の選択です'. The 'Which amplicon do you want to use in the instrument run?' section has a dropdown menu set to 'Fast (~ 40 minutes to complete a run)'. A callout points to this dropdown: '使用している試薬の添付文書で確認してください。'. A callout on the left side of the window points to the 'Experiment Properties' button in the 'Setup' menu: '“SYBR Green Reagent” を選択するとこのチェックボックスが出現します。'. The 'Run' button is visible in the top right corner.

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The screenshot shows the 'Define Targets and Samples' window in 7500 Software v2.0.6. The window title is 'Experiment: 20151214name etc Type: Standard Curve Reagents: TaqMan® Reagents'. The 'Define Targets' section has a table with columns: Target Name, Reporter, Quencher, and Colour. Two targets are listed: Target 1 and Target 2, both with Reporter FAM and Quencher NFQ-MGB. A callout points to this table: 'Target genes'. The 'Define Samples' section has a table with columns: Sample Name and Colour. Six samples are listed: Sample 1 through Sample 6, each with a different color. A callout points to this table: 'Samples'. The 'Assign Targets and Samples' button is visible at the bottom right. The 'Run' button is visible in the top right corner.

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